



# G-2030M Instruction

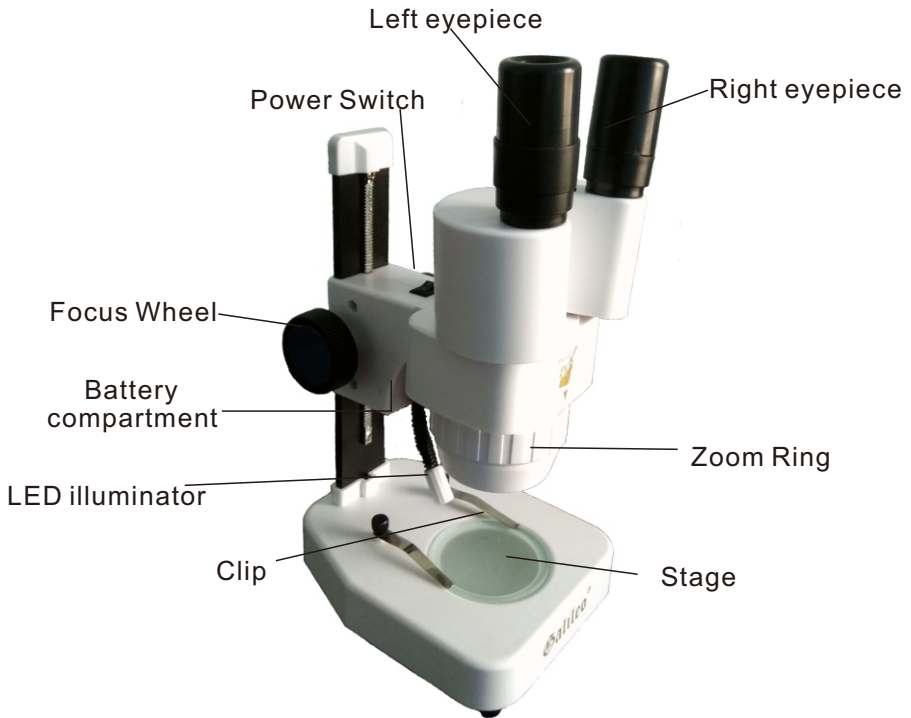

## How to use your stereomicroscope


1. Put the viewing object on the Stage, fixed with the clip.
2. Turn on the switch, the LED light is on.

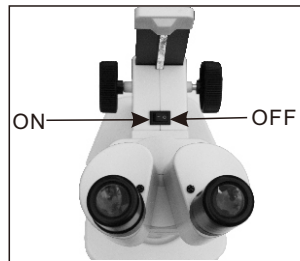


3. Adjust flexible LED light to proper position lighting object.



4. Set the interpupillary distance by grasping the eyepiece base and bending it until you can see one circle of view.

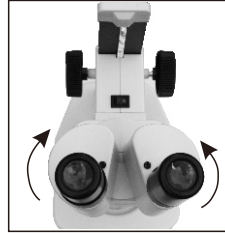

5. Close your left eye and sight the viewing object with your right eye, rotating the Focus wheel until the image is sharp and clear.

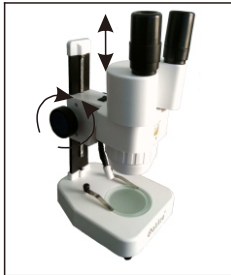

6. Open your left eye and rotating the left eyepiece until the viewing object is sharp and clear.




7. You can rotating the Zoom ring to change the magnification 20X or 30X.



8. Battery installation

Open the cover of Battery case, insert 2X AA batteries into the case, there is a indication +, - on the cover, then close the cover.

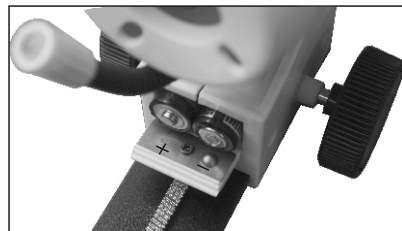

## SPECIFICATION

1. Magnification: 20X / 30X
3. Working distance: 80mm / 67mm
5. Illumination: LED
7. Weight: 780g

2. Field of view: 8mm / 5.5mm
4. Interpupillary distance: 52-74mm
6. Power: DC3V 2AA batteries